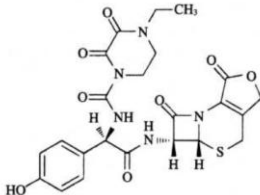
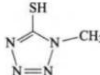
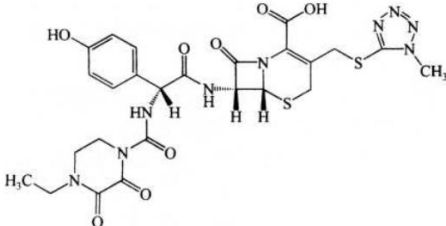
Test report

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Sample Information			
样品信息			
Sample Name 样品名称	cefoperazone 头孢哌酮	Sample properties 样品性状	White powder 白色粉末
Date of receipt 收样日期	2025/09/24	During testing 测试期间	2025/10/16~ 10/24
Test the ingredients and structural formula			
测试成分及结构式			
Cefoperazone, impurity A, impurity C, and isomer S. 头孢哌酮、杂质A、杂质C和异构体S. 杂质 A  (5aR, 6R) -6- [[(2R) -2- [[(4-乙基-2, 3-二氧代顺喹-1-基) 甲酰基] 氨基] -2- (4-羟基苯基) 乙酰基] 氨基] -5a, 6-二氢-3H, 7H-氮杂环丁二烯并 [2, 1-b] 咪唑并 [3, 4-d] [1, 3] 噻唑-1, 7 (4H) -二酮 杂质 c  1-甲基-5-巯基四氢唑 C2H4N4S 116.15 			
Experimental Requirements			
实验要求			
Method development. The separation between peaks in the system suitability chromatogram should be>1.5. The test sample solution should be injected continuously for 20 injections, and the retention time difference between the main peaks of the first and 20th injections should be <0.5 min. 方法开发。系统适用性色谱图中各峰之间的分离度应>1.5。供试品溶液连续进样 20 针，第 1 针与第 20 针主峰保留时间差应<0.5min。			

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Aqueous acetone solution (prepared by mixing 14 ml of triethylamine with 5.7 ml of glacial acetic acid, diluted to 100 ml with water, and shaken thoroughly) -acetonitrile-water mixture (1.2:120:880), with pH adjusted to 3.0 using glacial acetic acid.

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Solvent: Phosphate buffer: Take 39.0 ml of 0.2 mol/L sodium dihydrogen phosphate solution and 61.0 ml of 0.2 mol/L sodium dihydrogen phosphate solution, shake well, and adjust the pH to 7.0 with phosphoric acid.

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1.2.2. Test solution (供试品溶液):

Take cefoperazone about 25 mg, accurately weigh it, put it in a 50 ml volumetric flask, add a small amount of solvent to dissolve, then dilute with mobile phase to the mark, shake well, and get it.

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1.2.3. System suitability solution (系统适用性溶液):

Weigh approximately 2 mg of cefoperazone reference substance, impurity A reference substance (dissolved in acetonitrile first), impurity C reference substance, and cefoperazone S isomer into a 10 ml volumetric flask. Add a small amount of cosolvent to dissolve, then dilute with mobile phase to the mark and shake well to obtain the solution.

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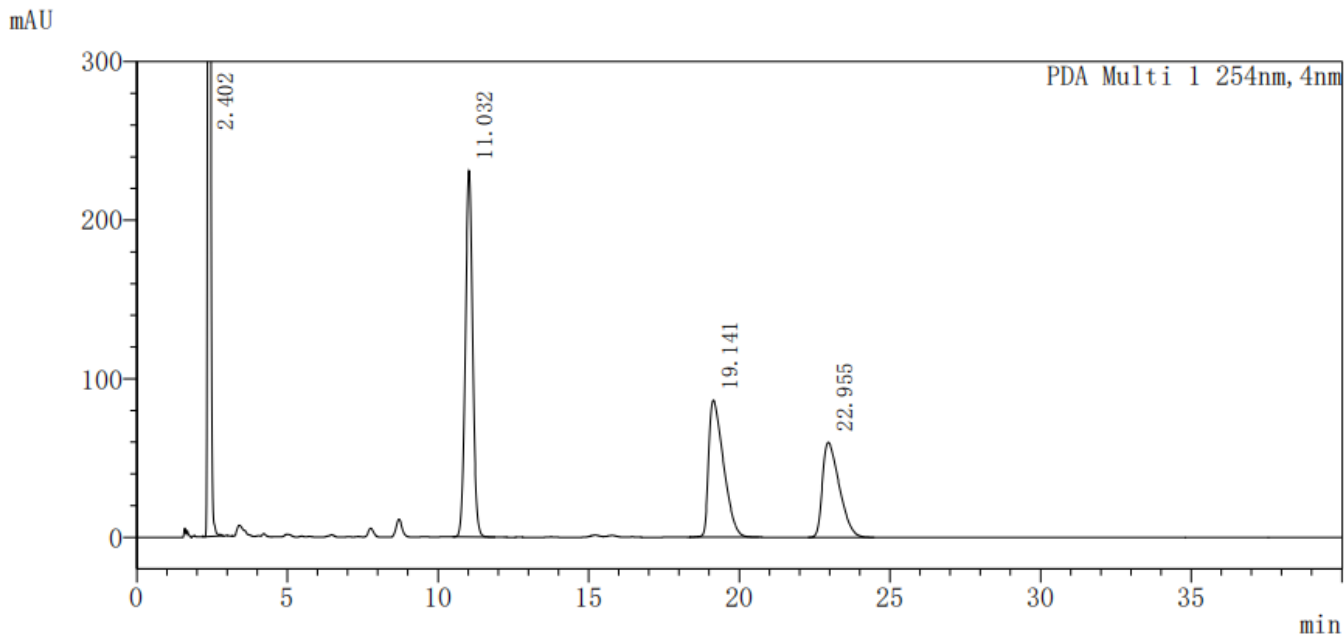
2. Chromatogram and data (谱图和数据)

(1) System suitability solution (系统溶液图)

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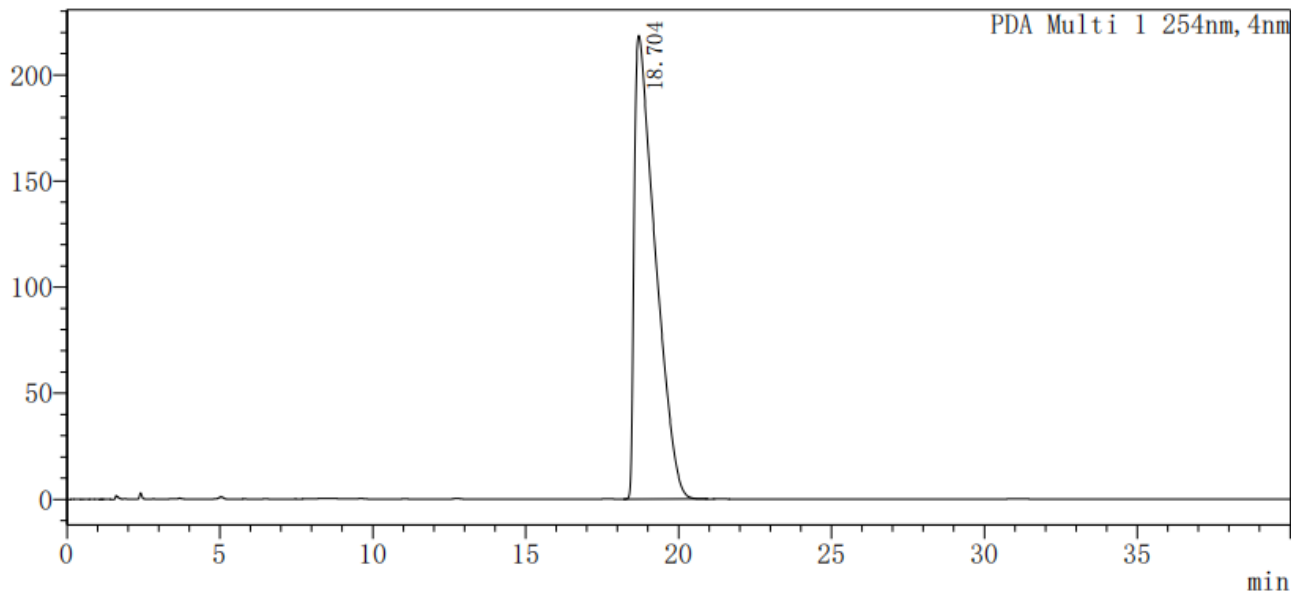
Conclusion (结论)

Using Ultimate[®] LP-C18 (4.6×150mm, 5μm, 300Å), the detection requirements can be met under these chromatographic conditions.

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