Test report

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Sample Information 样品信息			
Sample Name 样品名称	Ritagliptin 瑞他鲁肽	Sample properties 样品性状	White powder 白色粉末
Date of receipt 收样日期	2026/06/26	During testing 测试期间	2026/06/26~ 07/20
Test the ingredients and structural formula 测试成分及结构式			
H-Tyr-{Aib}-Gln-Gly-Thr-Phe-Thr-Ser-Asp-Tyr-Ser-Ile- $\{\alpha$ -Me-Leu}-Leu-Asp-Lys-{diacid-C20-gamma-Glu-(AEEA)-Lys}-Ala-Gln-{Aib}-Ala-Phe-Ile-Glu-Tyr-Leu-Leu-Glu-Gly-Gly-Pro-Ser-Ser-Gly-Ala-Pro-Pro-Pro-Ser-NH ₂			
Experimental Reuirements 实验要求			
Detection was performed using method kits and specialized columns, with amino acid residues falling within the specified range required by the client. 方法包和专用柱检测，氨基酸残基符合客户要求范围。			
Reference Method 参考方法			
方法包			
Reagent information 试剂信息			
Reagent Name 试剂名称	Grade 级别	Brand 品牌	
乙腈 产品编号: 00814-01010	HPLC	月旭	
乙酸	AR	国药	

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Report No (报告编号): 20260622-623

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1.1.1. Reference solution (对照溶液):

The two injections are designated as A and B respectively, diluted with a diluent to 1/5 of the original concentration. Take 160 μ L of the control solution from the method kit, transfer it to a test tube, and add 100 μ L of the diluted A solution and 100 μ L of the diluted B solution. Shake well and allow to react at room temperature for 60 minutes. Then add 400 μ L of n-hexane solution, tighten the cap, and shake for 5 - 10 seconds. Let stand at room temperature to allow layer separation. Collect 200 μ L of the lower layer, add 800 μ L of water, and mix thoroughly. Repeat the process with another 200 μ L of the lower layer and 800 μ L of water, mixing uniformly. Filter through a 0.22 μ m organic membrane before analysis.

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1.1.2. Test solution (供试品溶液):

Hydrolysis: Weigh precisely 50 mg of the test sample into a 10 mL volumetric flask, dissolve it in water, and dilute to the mark. Precisely transfer 1 mL of this solution to a 20 mL vial, add 9 mL of hydrolysis solution (prepared by accurately measuring 5.6 mL of hydrochloric acid into a 10.0 mL volumetric flask, diluting with water to the mark, and mixing thoroughly) to dissolve the sample; seal the vial under nitrogen atmosphere, place it in an oven at 110 $^{\circ}$ C for hydrolysis for 24 hours, remove it, allow it to cool to room temperature, then take 200 μ L of the solution, dry it under nitrogen, and dissolve it in 200 μ L of water.

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Derivation: The sample was diluted 50-fold with water. Absorb 160 μ L of the solution and transfer to a test tube. Add 100 μ L of diluted Solution A and 100 μ L of diluted Solution B, mix thoroughly, and incubate at room temperature for 60 minutes. Then add 400 μ L of n-hexane solution, tighten the cap, and shake for 5 - 10 seconds. Allow to stand at room temperature for stratification. Collect 200 μ L of the lower layer, add 800 μ L of water, and mix uniformly. Repeat the process with another 200 μ L of the lower layer and 800 μ L of water, mixing thoroughly. Filter through a 0.22 μ m organic membrane before analysis.

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